- Alkalinity Test Kit
- Trousse d'analyse alcalinité
- Alkalitäts-Test-Kit
- Kit para la determinación de alcalinidad

5–100, 20–400 mg/L

• Mod. AL-AP MG/L

• # 24443-01

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WARNING

Handling chemical samples, standards, and reagents can be dangerous. Review the Material Safety Data Sheets before handling any chemicals.

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Measuring Hints and General Test Information

- Wash all labware between tests. Contamination may alter test results. Clean with a non-abrasive detergent or a solvent such as isopropyl rubbing alcohol. Use a soft cloth for wiping or drying. Do not use paper towels or tissue on plastic tubes as this may scratch them. Rinse with clean water (preferably deionized water).
- When titrating, count each drop of titrant. Hold the dropper vertically. Swirl the mixing bottle after each drop is added.
- The result can be expressed in grains per gallon (gpg) by dividing the mg/L result by 17.1.
- To open PermaChem® Powder Pillows:
 1. Tap the bottom of the pillow on a hard surface.
 2. Tear open the pillow along the dashed line.
 3. Open the pillow and form a spout by squeezing the side edges.
 4. Pour the contents into the sample.
- Hach strongly recommends that, for optimum test results, reagent accuracy be checked with each new lot of reagents. Use the standard solution included in this kit or listed in the *OPTIONAL REAGENTS AND EQUIPMENT* section. Follow the instructions included with each standard solution.

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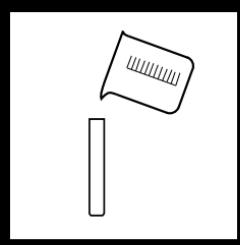
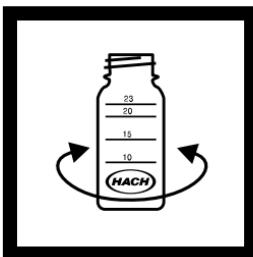
6. Multiply by 20 the number of drops of titrant used. This is the mg/L of phenolphthalein alkalinity as calcium carbonate (CaCO_3).

$\text{mg/L CaCO}_3 \text{ phenolphthalein alkalinity} = \text{number of drops} \times 20$

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7. Add the contents of one Bromcresol Green-Methyl Red Indicator Powder Pillow to the mixing bottle.

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8. Swirl to mix.

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- Agite para mezclar.

- 9.** Add Sulfuric Acid Standard Solution one drop at a time. Count each drop. Swirl the mixing bottle after each drop is added. Add drops until the sample turns pink.

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- 10.** Multiply by 20 the total number of drops of titrant used in both steps 5 and 9. This is the total mg/L of methyl orange alkalinity as calcium carbonate (CaCO_3).

$$\text{mg/L CaCO}_3 \text{ methyl orange alkalinity} = \text{number of drops} \times 20$$

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9. Multiply by 5 the total number of drops of titrant used in both steps 4 and 8. This is the total mg/L of methyl orange alkalinity as calcium carbonate (CaCO_3).

$\text{mg/L CaCO}_3 \text{ methyl orange alkalinity} = \text{number of drops} \times 5$

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REPLACEMENTS

Description	Unit	Cat. No.
Alkalinity Reagent Set,		
Drop Count Titration 0–400 mg/L as CaCO ₃	100 tests.....	24374-01
<i>Includes:</i> (1) 942-99, (1) 943-99, (1) 23497-32		
Bottle, mixing, glass.....	6/pkg.....	2327-06
Brom cresol Green-Methyl Red Powder Pillows.....	100/pkg.....	943-99
Instruction Card, AL-AP Test Kit	each.....	24443-89
Measuring Tube, plastic, 5.83 mL.....	each.....	438-00
Phenolphthalein Powder Pillows.....	100/pkg.....	942-99
Sulfuric Acid Standard Solution, 0.035 N.....	100 mL MDB ¹	23497-32

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